


Check the booking calendar. Are you the last user?

α 1

- 1•Remove your sample from the microscope
- 2•Clean the objective lens using lens cleaning tissue
- 3•Switch off the fluorescence bulb

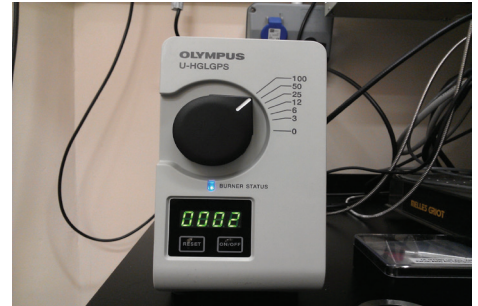

β 1

Turn off the Multiline argon (457/488/515) laser;

- 1•Turn the key anti-clockwise into the off position.
- 2•*WAIT* for 5 mins.
- 3•When the air blowing from the fan is cool
- 4•Turn the power SWITCH off.

WAIT

β 2




γ 1

Turn off the 559 laser;

- 1•Turn the key anti-clockwise to the off position.
- 2•Turn the power SWITCH off
(there's no need to wait here!)

γ 2

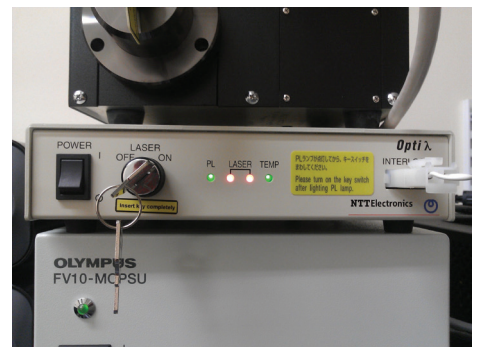

δ 1

Turn off the microscope keypad.
The On/Off button is on the BACK.



ϵ 1

Log off the computer.

ϵ 2

Switch off the controller box
(the black button on the left)

