

Live Cell Olympus Turn ON

A1

Switch on the controller boxes
(the black buttons on the left)

A2



B1

Turn on the 559 laser;

1•turn the power SWITCH on.

2•The TEMP LED (green) will flash.

3•*WAIT*

4•After approximately 2 mins the LED will stop flashing.

5•Only then turn the KEY clockwise to the on position.

6•The LASER LEDs (red) will flash

7• *WAIT*.

8•When these red LEDs stop flashing the laser is ready to use

WAIT

B2



C1

Turn on the Multiline argon (457/488/515) laser;

1•turn the power SWITCH on.

2•Then turn the key clockwise to the on position
(there is no need to wait for anything here!).

C2



D

Turn on the fluorescence bulb



E

Turn on the microscope keypad.
The On/Off button is on the BACK.

It's crucial to do this before starting the software

