

# Olympus SoRa Turn OFF



**Check the booking calendar. Are you the last user?**

**$\alpha 1$**

- 1•Remove your sample from the microscope
- 2•Clean the objective lens using lens cleaning tissue
- 3•Close software
- 4.Shut down PC

**$\beta$**

Turn off the microscope keypad.  
The On/Off button is on the BACK.



**$\gamma$**


Turn off the spinning disc;



**$\delta$**

Turn off the lasers;

- 1•Turn the key anti-clockwise into standby position.
- 2•Turn the power SWITCH off.

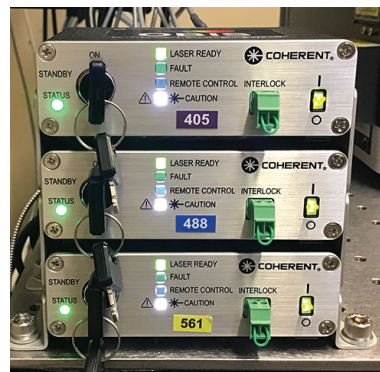

**$\epsilon 1$**

**$\epsilon 2$**

**$\epsilon 3$**

Switch off the controller box  
(the orange buttons on the left)

