

A3

The image shows three Coherent laser units stacked vertically. Each unit has a 'STANDBY' label with a green indicator light that is illuminated. To the right of the 'STANDBY' label is a 'STATUS' label with a green indicator light that is also illuminated. Each unit has a 'LASER READY' label with a green indicator light that is illuminated. To the right of the 'LASER READY' label is a 'FAULT' label with a red indicator light that is not illuminated. Each unit has a 'REMOTE CONTROL' label with a green indicator light that is illuminated. To the right of the 'REMOTE CONTROL' label is a 'CAUTION' label with a yellow indicator light that is illuminated. Each unit has a 'COHERENT' logo. The units are labeled with their wavelengths: 405, 488, and 561. Each unit has a green safety interlock connector. The units are connected to a power supply and a ground connection.

C2

E

Turn on the microscope keypad.
The On/Off button is on the BACK.
It's crucial to do this before starting the software.