



A. Switch on the two controller boxes
(the black buttons on the left)

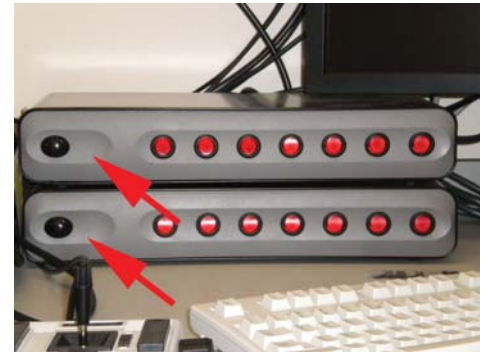


B. Turn on the 559 laser;
1•turn the power SWITCH on.
2•The TEMP LED (green) will flash.
3•*WAIT*
4•After approximately 2 mins the LED will stop flashing.
5•Only then turn the KEY clockwise to the on position.
6•The LASER LEDs (red) will flash
7• *WAIT*.
8•When these red LEDs stop flashing the laser is ready to use



C. Turn on the Multiline argon (457/488/515) laser;
1•turn the power SWITCH on.
2•Then turn the key clockwise to the on position
(there is no need to wait for anything here!).

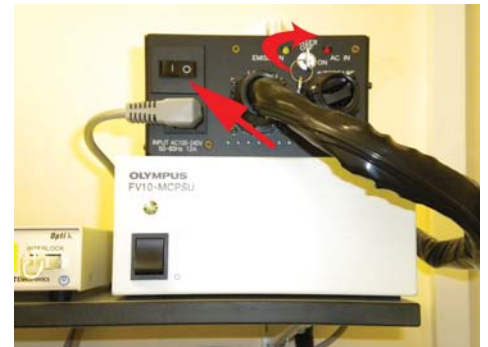

D. Turn on the fluorescence bulb

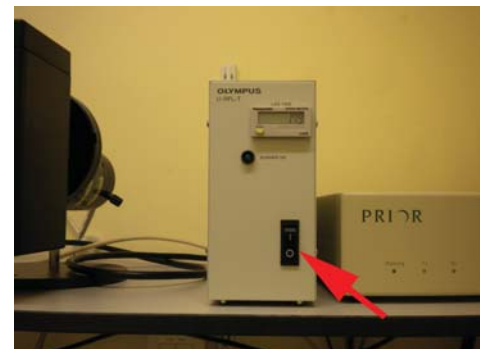

E. Start the program;
1•double click the icon 'FV10-ASW'

